## IKA® RW 47 D

Indicaciones de seguridad	ES	27
Veiligheidsinstructies	NL	29
Norme di sicurezza	IT	30
Säkerhetsanvisningar	SV	32
Sikkerhedshenvisninger	DA	34
Sikkerhedsanvisninger	NO	35
Turvallisuusohjeet	FI	37
Normas de segurança	PT	38
Υποδείξεις ασφαλείας	EL	40

Betriebsanleitung	DE	3
Operating instructions	EN	11
Mode d'emploi	FR	19

Wskazówki bezpieczeństwa	PL	42
Bezpečnostní upozornění	CS	44
Biztonsági utasítások	HU	45
Varnostna navodila	SL	47
Bezpečnostné pokyny	SK	49
Ohutusjuhised	ET	50
Drošības Norādes	LV	52
Saugos Reikalavimai	LT	54
Инструкции за безопасност	BG	55
Indicații de siguranță	RO	57

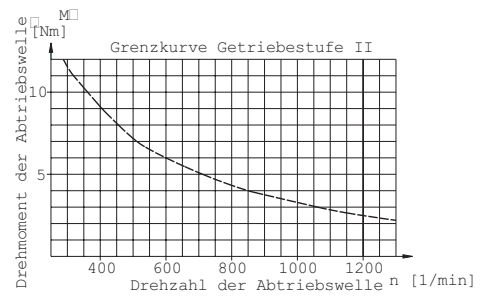
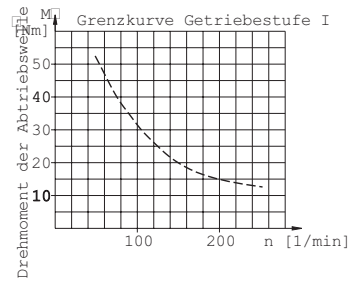
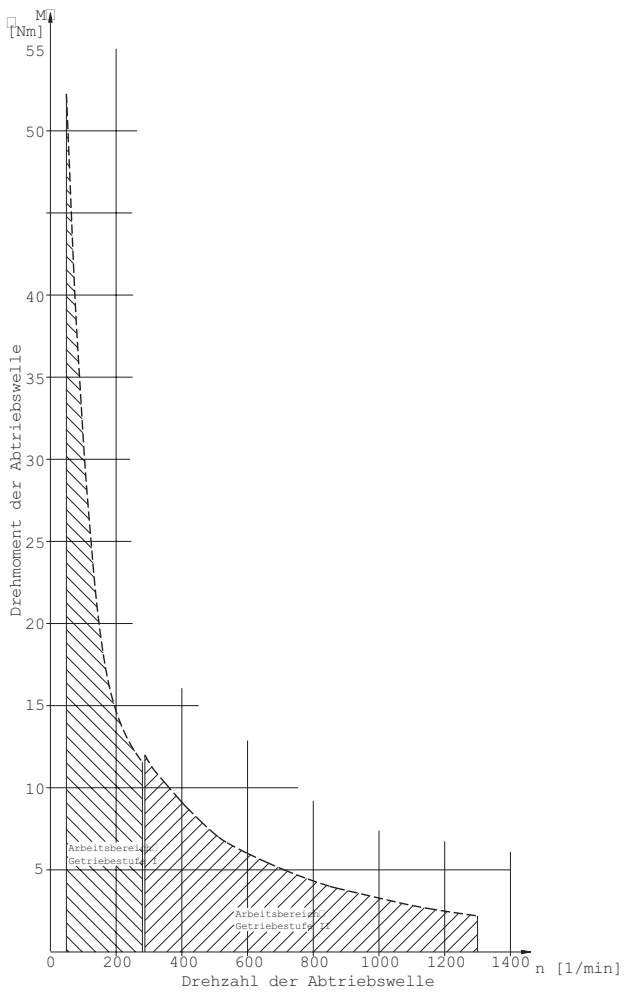
Fig. 1

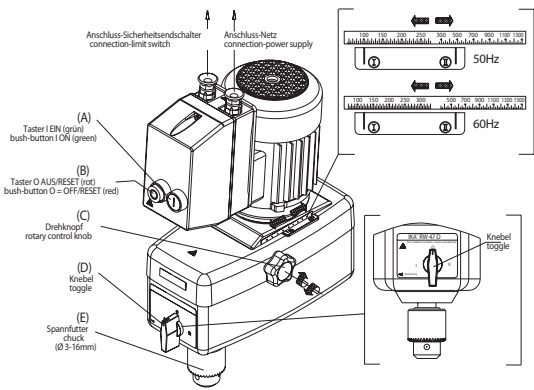
Fig. 2

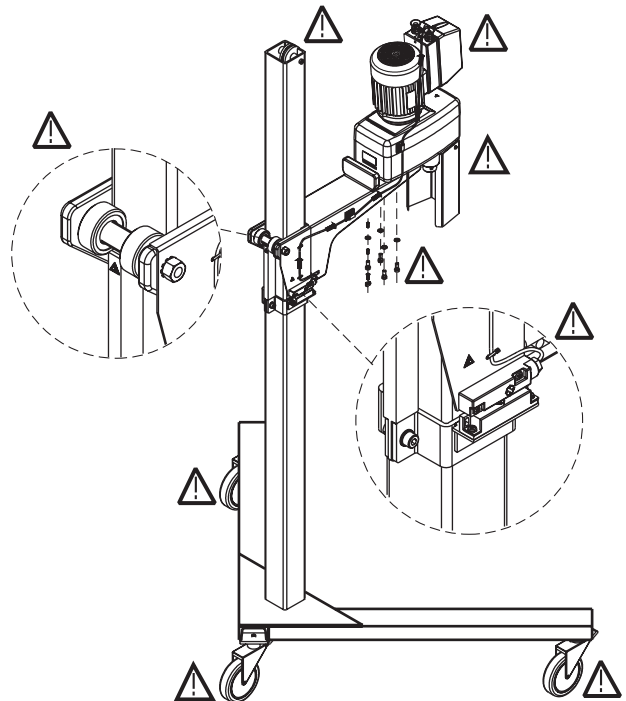
Fig. 3

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## Contents

	Page
CE-Declaration of conformity	11
Explication of warning symbols	11
Safety instructions	12
Unpacking	13
Correct use	13
Useful information	14
Drive	14
Motor circuit breaker	14
Speed display	15
Gear speed change	15
Commissioning	15
Switching on the device	15
Output shaft	16
Maintenance and cleaning	16
Permitted <b>IKA</b> <sup>®</sup> stirring tools	16
Accessories	17
Technical data	17
Troubleshooting	18
Warranty	18
Wiring diagram	59

## CE - Declaration of conformity

We declare under our sole responsibility that this product corresponds to the regulations 2006/42/EEC and 2004/108/EEC and conforms with the standards or standardized documents: DIN EN IEC 61 010-1, -2-051; DIN EN ISO 12 100-1, -2; DIN EN IEC 61 326-1 and EN 60 204-1.

## Explication of warning symbols

General hazard

This symbol identifies information **that is of vital importance for safeguarding your health and safety**. Disregarding this information can lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the appliance**. Disregarding this information can result in damage to the appliance.

This symbol indicates information **which is important for ensuring that the operations of the appliance are performed efficiently and for using the appliance**. Failure to observe this information can result in inaccurate results.

## Safety instructions

### *For your protection*

- **Read the operating instructions in full before starting up and follow the safety instructions.**

- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise there is a risk of:

- splashing liquids
- projectile parts
- body parts, hair, clothing and jewellery getting caught.

- Uncontrolled reactions can be triggered by the energy generated. In case of increased operational hazards, users must take additional appropriate safety precautions (e.g. shatter protection). In any case, **IKA®** recommends that users who use critical or hazardous materials in their processes use additional appropriate measures to ensure safety in the trial set-up. For example, users can implement measures that inhibit fire or explosions or comprehensive monitoring equipment.
- The device must be wired by a qualified technician before first use in order to prevent risk of electric shock!

- Please pay attention to the dangerous parts of the equipment, see fig. 3.
- Use a stirring shaft protective equipment and a safety cutout for height adjustable stands!
- The high torque developed by the **IKA®** RW 47 D requires particular care in the choice of stand, cross sleeve and anti-rotation element for the agitating vessel.
- If the emergency stop switch is not within reach when the unit is operating, an additional emergency stop switch must be installed within easy reach, in the working area.
- Before commissioning, the correct direction of rotation of the motor must always be ensured (test run without stirring tool: motor rotation as indicated on the arrow plate, or clockwise rotation, looking down on motor). The wrong direction of rotation can lead to the stirring tool coming off the shaft.

- If the unit is operated in different locations with a 5-pin plug, the direction of rotation must be checked before commissioning, with no stirring tool fitted.
- The appliance is not suitable for manual operation.
- Set up the stand in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface
- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
- The agitated vessels used for stirring have to be secured. Consider on a good stability of the entire structure.
- Secure the stirring vessel against twisting.
- Never operate the equipment with the stirrer tools rotating freely.
- You must ensure that the stirring element is securely clamped in the chuck!
- The stirring elements may only be inserted into the hollow shaft to such a depth that they do not abrade the top edge of the casing.
- Please observe the permitted speed for the stirring element. Never set higher speeds.
- Make certain that the unit is set at the lowest speed before commissioning; otherwise, the unit will begin running at the speed last set. Gradually increase the speed.
- Pay attention when setting the speed to any imbalance of the stirrer tools and possible spraying of the medium to be stirred.
- Reduce the speed if
  - the medium splashes out of the vessel because the speed is too high
  - the appliance is not running smoothly
  - the appliance begins to move around because of dynamic forces
- The unit must be switched off (motor standstill) before changing to another gear speed, otherwise the toothed gears in the reduction stage will be damaged.
- Do not touch rotating parts during operation!!
- Avoid knocks and impacts on the lower end of the shaft and the chuck gear teeth. Even minor, invisible damage can lead to imbalance and uneven shaft action.
- Ensure that the stand does not start to move.
- It is important to note that the surfaces of the motor (cooling fins) and certain parts of the bearing can get very warm during operation.

- Never cover the ventilation slots or cooling fins on the motor as these are needed for cooling the drive system.
- Imbalance of the output shaft, the chuck and in particular the stirring tools can lead to uncontrolled resonant vibrational behaviour of the equipment and the whole assembly. Glass apparatus and stirrer containers can be damaged or shattered by this. It can cause injury to the operator, as can the rotating stirring tool. In this case exchange the stirring tool for one without imbalance or remedy the cause of the imbalance. If there still is imbalance, return it to the dealer or the manufacturer along with a description of the fault.
- Beware of the risk of
  - flammable materials
  - glass breakage as a result of mechanical shaking power.
- There may be electrostatic activity between the medium and the output shaft which could pose a direct danger.
- Process pathogenic materials only in closed vessels under a suitable extractor hood. Please contact **IKA®** if you have any questions.
- Do not operate the appliance in explosive atmospheres, with hazardous substances or under water.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- Safe operation is only guaranteed with the accessories described in the "Accessories" chapter.
- Always disconnect the plug before fitting accessories.
- The appliance can only be disconnected from the mains supply by pulling out the mains plug or the connector plug.
- The socket for the mains cord must be easily accessible.
- After an interruption in the power supply, or a mechanical interruption during a stirring process, the unit does not automatically restart.

#### *For protection of the equipment*

- The appliance may only be opened by experts.
- The voltage stated on the type plate must correspond to the mains voltage.
- Socket must be earthed (protective ground contact).
- Protect the appliance and accessories from bumps and impacts.

## Unpacking

- **Unpacking**
  - Please unpack the device carefully
  - In the case of any damage a detailed report must be set immediately (post, rail or forwarder)
- **Delivery scope**  
**Stirring unit IKA® RW 47 D**
  - IKA® RW 47 D stirring unit as per order,
  - Chuck key,
  - Allen key,
  - 4 Allen screws
  - Operating instructions

## Correct use

- **Use**  
The laboratory stirrer **IKA®** RW 47 D can be used to stir and mix liquids of low to medium viscosity with various stirring tools.
- Operating modes: on stand

### • Range of use

- Laboratories
- Pharmacies
- Technical colleges
- Small-scale production

The safety of the user cannot be guaranteed if the appliance is operated with accessories that are

not supplied or recommended by the manufacturer or if the appliance is operated improperly contrary to the manufacturer's specifications or if the appliance or the printed circuit board are modified by third parties.

## Useful information

You can set speeds of 57 to 1300 rpm with the 50 Hz (mains frequency) version and speeds from 69 to 1560 rpm with the 60 Hz (mains frequency) version. The unit heats up during use. The ample cooling surfaces on the motor serve to distribute and radiate the heat as evenly as possible.

You can also use high-sided mixing vessels together with **IKA**® stands without any problem. The unit comes without a connection cable because the user's installations mean that it is not

possible to issue a standard length cable due to the different sizes of premises. Experts can determine the pin assignment from the wiring plan.

The precision shaft is shaped as a hollow shaft and accommodates stirrer shanks of up to 13 mm in diameter. It is important to note here that these may only be inserted to just before the top edge of the casing. You must also ensure that the stirrer shank is securely clamped in the chuck.

## Drive

The motor can be infinitely adjusted via a friction wheel drive unit but is operated in one position only. The power output, rotational speed and torque of the motor are regarded as constant and are optimized for this operating position. After the friction wheel drive the motor power output is transmitted to the off-drive shaft of the stirrer via either the first or second stage of the subsequent manually adjustable toothed gear train.

Neglecting all losses, the power at the off-drive shaft during power transmission (conversion) is always as great as the power output of the

motor. The gear train merely provides rpm and torque conversion, which can be described in purely theoretical terms by the characteristic curve plotted in Fig. 1.

Friction wheel wear is kept low by a helical clutch matched to the gear train. The down-pressure required at the friction wheel is adjusted by the helical clutch according to the moment acting on the stirrer shaft. A low torque results in a low down-pressure; a high torque results in a high downpressure.

## Motor circuit breaker

The ball race three-phase motor is maintenance-free. Its attached switching unit features a special safety device for operation in the case of over-current and undervoltage (undervoltage release) to permanently disconnect the motor and avoid thermal damage. The appliance can only be started up again by activating the ON/OFF switch (Fig. 2), when power permitted for the motor is supplied.

You can find the basic connection and rating for the motor on its rating plate.

The safety switch SI 400 can also be connected in the motor switching unit. When used in combination with an **IKA**® stand, this allows the user to switch on the appliance only at a defined height in the stand's working range.

If the stirring unit is operated in such a way that the appliance's EMERGENCY OFF switch ( (B) (Fig. 2) is not accessible, you must add another EMERGENCY OFF switch to the work area so that it can be easily reached.

## Speed display

The speed on the **IKA**® RW 47 D agitator is adjusted by means of the rotary button (C). You can read the set speed from the marking on

the top of the casing on the label attached to the motor support. (The values are only valid for low agitator shaft lows).

## Gear speed change

You can select one of the three gear positions available (0, I or II) using the toggle button (D) on the front of the appliance. You may only switch to a different cog wheel gear

level when the appliance is switched off (motor stopped). Otherwise this may cause damage to the gear cogs. The arrow on the front plate indicates the direction of the chuck.

## Commissioning

The laboratory stirrer **IKA**® RW 47 D must be secured to a stable stand with four screws (e.g. R 474 or R 472 ) and wired according to the wiring plan for proper use. The stirring bowl must always be well secured for safety reasons. You must also ensure that the mounting equipment (e.g. floor stand, telescope stand) is set up in such a way that it is not liable to topple and does not start to move during the stirring procedure. Accessories must be assembled and wired according to the assembly instructions.

If the stirring unit is operated in such a way that the appliance's EMERGENCY OFF switch (B) (Fig. 2) is not accessible, you must add another EMERGENCY OFF switch to the work area so that it can be easily reached. Points of particular danger (such as crushing, rotating parts, clamping screws for stand legs and rollers etc.) are marked in Fig. 3 with an exclamation mark. Please pay particular attention to these danger points during operation!

## Switching on the device

Check whether the voltage given on the type plate corresponds to the available mains voltage. The socket used must be earthed (fitted with earth contact). If these conditions have been met, the machine is ready to operate when the mains plug is plugged in. If these conditions are not met, safe operation is not guaranteed and the machine could be damaged.

When working with safety switch SI 400, the appliance only starts up once it has been set to the designated working height and the switch contact activates the safety switch.

During commissioning of the machine the output shaft starts to run at the last speed set. Therefore check the setting of the control knob (speed).

Also ensure that the speed set is suitable for the test texture selected. If you are unsure, set the motor slide to the lowest speed front motor slide stop) using the side knob and the gear to the lowest level (Fig. 2).

The appliance will start up when you press the ON switch on the motor (Fig. 2).

If the unit has not been used for some time, a knocking noise will be heard when switching on, which is due to the preload on the friction lining of the friction wheel drive. This has no detrimental effect on the function of the unit, and the knocking noise will disappear after a short running-in period.

## Output shaft

The chuck (E) (Fig. 2) and off-drive shaft allow stirring elements with shank diameters of 3 mm to 16 mm to be clamped. Shank diameters of up to 13 mm can be inserted through the hollow shaft

to just before the top edge of the inside casing. You must ensure that the stirring shank is securely clamped in the chuck.

## Maintenance and cleaning

The stirrer is maintenance-free.

Only clean **IKA®** appliances using these **IKA®** approved cleaning agents:

Dirt	Cleaning agent
Dyes	Isopropanol
Building materials	Water containing detergent, Isopropanol
Cosmetics	Water containing detergent, Isopropanol
Food	Water containing detergent,
Fuels	Water containing detergent,
Other materials	Please consult <b>IKA®</b>

- Wear protective gloves when cleaning the devices.
- Do not place electrical appliances into the cleaning agents for cleaning purposes.
- Do not allow moisture to get into the appliance when cleaning.
- Please consult **IKA®** before using any cleaning or decontamination methods, other than those recommended here.

- When cleaning the unit, take particular care not to bring substances containing oil or grease into contact with the running surfaces of the cone pulley and the friction wheel. This would substantially reduce the frictional coefficient resulting from the pairing of the friction wheel and cone pulley materials, and would impair power transmission in the unit.
- Make certain that the hexagon head nut (part no. 153 of spare parts diagram) is sufficiently tightened in order to prevent the shifting of the motor plate and thus the speed as well.

### Spare parts order

When ordering spare parts, please give:

- Machine type
- Manufacturing number, see type plate
- Item and designation of the spare part, see **www.ika.com**, spare parts diagram and spare parts list

### Repair

In case of repair the device has to be cleaned and free from any materials which may constitute a health hazard.

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

## Permitted IKA® Stirring tools

Max. speed rpm

R 2305	Propeller stirrer 4fl	1300	R 1302	Dissolver stirrer	1000
R 2302	Propeller stirrer 4fl	600	R 1303	Dissolver stirrer	2000
R 1385	Propeller stirrer 3fl	800	R 1376	Surface stirrer	800
R 1388	Propeller stirrer 3fl	400	R 2311	Surface stirrer	600
R 1313	Turbine stirrer	800	R 1333	Anchor stirrer	800
R 1300	Dissolver stirrer	2000	R 1393	Worm stirrer	1000
R 1301	Dissolver stirrer	800			

## Accessories

R 472	Floor stand	SI 400	Safety limit switch
R 474	Telescopic stand	SI 472	Support holder
R 302	Stirring shaft protection	SI 474	Support holder

## Technical data

### Speed range

50 Hz stage I	rpm	57-275
50 Hz stage II	rpm	275 – 1300
60 Hz stage I	rpm	69 - 330
60 Hz stage II	rpm	330 – 1560
Speed adjustment		Mechanical rotary button on the side
Speed display		Speed label on the side of the device
Max. torque stirrer shaft (overload)	Ncm	4642 ( bei 57 rpm )
Max. torque stirrer shaft	Ncm	3000 ( bei 100 rpm )
Max. torque stirrer shaft	Ncm	285 ( bei 1000 rpm )

### Motor data

Design voltage	VAC	3 x 400 Y	3 x 230 Δ	3 x 400 Y	3 x 230 Δ
Design frequency	Hz	50	50	60	60
Motor speed	rpm	2880	2850	3425	3425
Input power	W	554	548	489	490
Output power	W	370	370	370	370
Maximum output at stirrer shaft	W	300			
Permitted on-time	%	100			
Protection type to DIN EN 60529		IP 44			
Oversvoltage category		II			
Protection class		I (Protective earth)			
Contamination level		2			
Protection on overload		overcurrent and undervoltage switch on the three-phase motor			
Perm. ambient temperature	°C	+ 5 bis + 40			
Perm. humidity (rel.)	%	80			
Operating position		On stand, clamping chuck pointing down			
Drive		Rib cooled three-phase motor with friction wheel drive and subsequent 2-stage toothed gear train			
Maximum stirring quantity (water)	l	200			
For high-viscosity (HV)	mPas	10000-100000			
Clamping chuck - clamping range	mm	3-16			
Hollow shaft internal diameter	mm	13 (upto short before the internal housing upper edge)			
Stand attachment		Flange on housing lower part			
Housing material		Al, hermoplastic			
Dimension without extension arm (WxDxH)	mm	145x445x340			
Weight	kg	15			
Operation at a terrestrial altitude	m	max. 2000			

*Subject to technical changes!*

## Troubleshooting

In cases of fault or overload operation the device can become very hot (motor surface, cooling fins, bearings).

If the device is operated too long in overload or if the ambient temperature is too high, the device switches off permanently.

In case of fault, proceed as follows:

- Switch the device off (power supply OFF) and allow the device to cool down.
- Remove the stirrer tool and remove the device from the assembly.
- Reduce the speed and switch on the device without the stirrer tool (power supply ON).

If the actions described fail to resolve the fault

- Contact the service department,
- Send the device for repair, including a short description of the fault.

## Warranty

In accordance with **IKA**<sup>®</sup> warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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IKA® India Private Limited
IKA® Works Inc.