



MC 905 Laboratory detergent

Concentrated detergent and disinfecting solution for thorough cleaning of metallic, plastic and glass instruments, reservoirs, bottles and parts. Recommended for use in doctor practices and laboratories of any type.

Preparation / application:

- Prepare a 1-3% solution (10 - 30 mL / L)
- Use a 5% solution for highly soiled items
- Immerge items to be cleaned completely in the solution
- Leave at room temperature for at least 30 minutes
- Thoroughly rinse with distilled water
-

Sterilize cleaned items when needed

Ordering information

Cat. No.

MC001

Description

MC 905 liquid

Packaging

950 mL